


Installation Guidelines


LABORATORY/UTILITY SINKS

Small Model C5201
Large Model C5241



General Dimensions

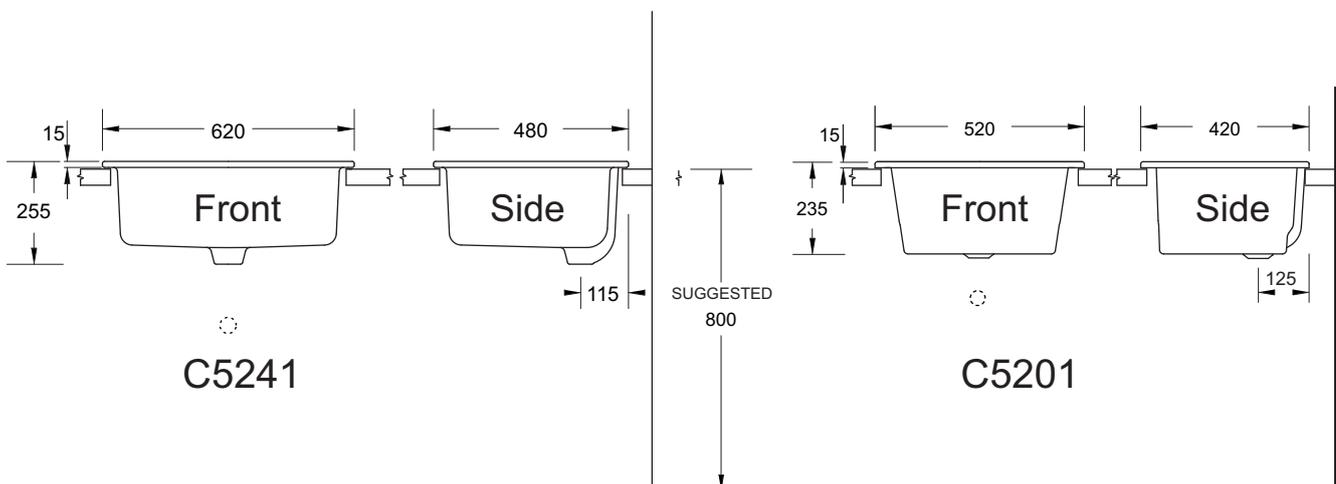


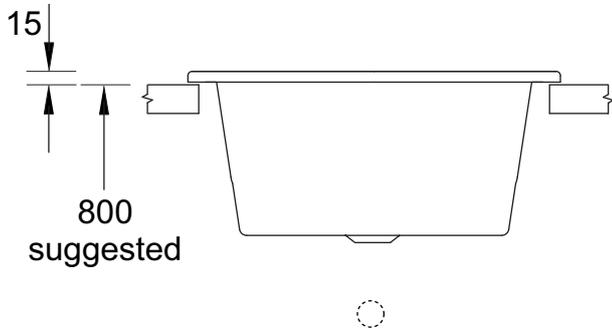

INSTALLATION


Inset Installation

- Place the sink at the required position.
- Mark the cutout in the benchtop in that position and according to supplied template.
- Cut benchtop.
- Place the sink in place, and seal around the sides with a suitable silicone.

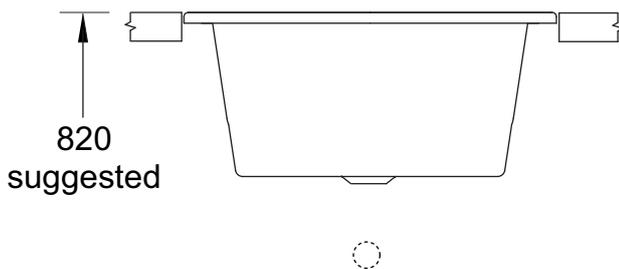

Flush Installation

- Place the sink at the required position.
- Mark the cutout in the benchtop using the sink as a template.
- Cut benchtop.
- Place the sink in place, and support from the underside.
- Seal around the sides with a suitable silicone.

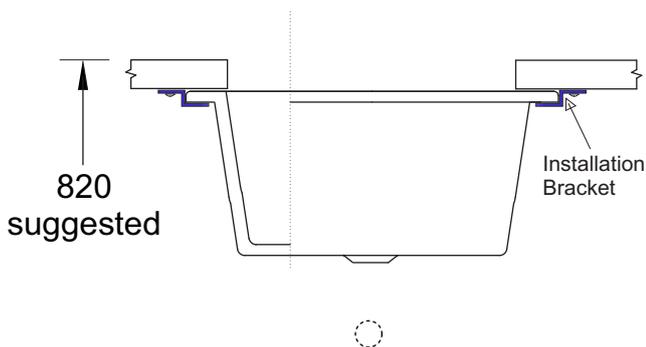

Undermount Installation

- Place the sink at the required position.
- Mark the cutout in the benchtop using the internal dimensions of the sink.
- Cut benchtop.
- Apply construction grade adhesive to the top rim of the sink, secure the sink to the underside of the benchtop, and allow to cure.
- add 4 support brackets on the underside.
- Seal around the inside with a suitable silicone.